



Alcohol Testing Program

INSTRUMENT PROCESSING SHEET

Agency Quincy PD S/N 80-000965
Date In 7/27/17 Date Out 8/18/2017
Ship P/U H/D CMI EE

Intake Performed By [Signature]
Registration
Annual
Return from CMI
Return from Enforcement Electronics
Other
Visual Inspection:
Case Handle
Dry Gas Holder Feet
Keyboard/Plug Back/Plugs
Screws tight Breath Hose
Other Equipment:
Power cord
Printer Cable
Other

Quality Checks Performed By [Signature]
Lab Temp 22.5
Breath Tube Screen
Replace External O-Rings
Instrument Set Up Verified
R-Value 176
Flow Verification (L/s)
Flow Column # ATP102
32mm 1160 (.139 - .169)
36mm 1171 (.156 - .190)
53mm 1246 (.228 - .278)
103mm 1498 (.447 - .547)
Barometric Pressure Check
Gauge ID # 28421
Stability Checks
Table with Simulator, Serial #, Lot #/Exp

Flow Calibration Performed By [Signature]
Flow Calibration N/A
Flow Calibration Complete
Flow Column #
5L/min 47mm
15L/min 53mm
30L/min 103mm
R-Value
Post Calibration Verification (L/s)
Flow Column #
32mm (.139 - .169)
36mm (.156 - .190)
53mm (.228 - .278)
103mm (.447 - .547)
Maintenance Performed By
Battery Replacement
Dry Gas Regulator Replacement
Breath Tube Replacement
Other
Quality Checks Cont. Performed By [Signature]
Simulator Temperatures 34C
External Digital Therm. ID# 300505
34C +/- .2 Serial #: SD3962
34C +/- .2 Serial #: SD1032
34C +/- .2 Serial #: SD3933

RECEIVED
AUG 21 2017
FDLE
Alcohol Testing Program

Calibration Adjustment Performed By
Calibration Adjustment N/A
Calibration Adjustment Complete
Barometric Pressure Gauge ID #
Table with Simulator, Serial Number, Lot Number, Expiration
Post Calibration Adjustment Stability Checks
Table with Simulator, Serial Number, Lot Number, Expiration

Department Inspection Performed By [Signature]
Barometric Pressure 1013 Gauge
ID# 28421 1013 Instrument
Mouth Alcohol Solution Lot # 2016-C
Acetone Stock Solution Lot # 2017-A
Table with Simulator, Serial Number
Attachments
Form 41
Pre-Stability Tests
Flow Calibration
Calibration Adjustment
Post-Stability Tests
Other

Notes/Suggested Service:
PA/AC OK SP
[Signature]
Date 8/21/17

Instrument Complies with Chapter 11D-8, FAC
Instrument Does Not Comply with Chapter 11D-8, FAC
Return to/Place into Evidentiary Use
Remain Out of Evidentiary Use
Conduct an Agency Inspection Before Evidentiary Use

Quality Control Review

QUINCY PD
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000965
08/18/2017
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	10:28
Control Test	0.049	10:28
Air Blank	0.000	10:29
Control Test	0.048	10:30
Air Blank	0.000	10:30
Control Test	0.048	10:31
Air Blank	0.000	10:31
Control Test Stats		
Average	0.0483	
Std Dev	0.0006	
Rel Std Dev(%)	1.1945	

P Murphy

Operator's Signature

QUINCY PD
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000965
08/18/2017
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	10:22
Control Test	0.079	10:23
Air Blank	0.000	10:24
Control Test	0.080	10:24
Air Blank	0.000	10:25
Control Test	0.080	10:25
Air Blank	0.000	10:26
Control Test Stats		
Average	0.0797	
Std Dev	0.0006	
Rel Std Dev(%)	0.7247	

P Murphy

Operator's Signature

QUINCY PD
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000965
08/18/2017
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	10:34
Control Test	0.198	10:35
Air Blank	0.000	10:35
Control Test	0.197	10:36
Air Blank	0.000	10:36
Control Test	0.199	10:37
Air Blank	0.000	10:37
Control Test Stats		
Average	0.1980	
Std Dev	0.0010	
Rel Std Dev(%)	0.5051	

P Murphy

Operator's Signature

QUINCY PD
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000965
08/18/2017
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	10:48
Control Test	0.080	10:48
Air Blank	0.000	10:49
Control Test	0.079	10:49
Air Blank	0.000	10:50
Control Test	0.080	10:50
Air Blank	0.000	10:50
Control Test Stats		
Average	0.0797	
Std Dev	0.0006	
Rel Std Dev(%)	0.7247	

DGS

P Murphy

Operator's Signature

SP

8/21/17
SP