



INSTRUMENT PROCESSING SHEET

Agency Madison County SOS/N 80-000779Florida Department of
Law EnforcementDate In 5/13/25DI Completion Date 5/30/2025☐ Ship ☒ P/U ☐ H/D ☐ CMI ☐ EE

Intake	By SP	Date <u>5/20/25</u>	Quality Checks	By PN	Date <u>5/30/2025</u>	Flow Calibration	By	Date																																																																	
☒ Annual ☐ Registration ☐ Return from CMI / EE Visual Inspection: ☒ Case ☐ Handle ☒ Keyboard ☒ Dry Gas Shelf ☒ Feet ☐ Breath Tube ☒ Ports ☐ Screws Tight Other Equipment/ Accessories: ☐ Power cord ☐ Printer Cable ☐ Static Bag ☐ 12V DC Cable Notes: AI indicates instrument needs battery change.			☒ Breath Tube Screen ☒ Replace External O-Rings ☒ Instrument Set Up Verified ☒ R-Value <u>230</u> ☒ Flow Verification (L/s) Flow Column # <u>ATP 105</u> 32 mm <u>.156</u> (.139 - .169) 36 mm <u>.171</u> (.156 - .190) 53 mm <u>.242</u> (.228 - .278) 103 mm <u>.527</u> (.447 - .547) ☒ Barometric Pressure Check Gauge ID # <u>28427</u> ☒ Stability Checks			Flow Column # _____ ☐ 5L/min – 17mm ☐ 15L/min – 53mm ☐ 30L/min – 103mm ☐ R-Value _____ ☐ Post Calibration Verification (L/s) Flow Column # _____ 32 mm _____ (.139 - .169) 36 mm _____ (.156 - .190) 53 mm _____ (.228 - .278) 103 mm _____ (.447 - .547)																																																																			
			<table border="1"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #/Exp</th></tr></thead><tbody><tr><td>0.050</td><td>MP5088</td><td>202406K 06/19/2026</td></tr><tr><td>0.080</td><td>MP5089</td><td>202406L 06/19/2026</td></tr><tr><td>0.200</td><td>MP5090</td><td>202406N 06/20/2026</td></tr><tr><td>0.080 DGS</td><td>N/A</td><td>AG429602 10/22/2026</td></tr></tbody></table>			Simulator	Serial #	Lot #/Exp	0.050	MP5088	202406K 06/19/2026	0.080	MP5089	202406L 06/19/2026	0.200	MP5090	202406N 06/20/2026	0.080 DGS	N/A	AG429602 10/22/2026	<table border="1"><thead><tr><th colspan="2">Maintenance By PN</th><th>Date <u>5/29/2025</u></th></tr></thead><tbody><tr><td colspan="2">☒ Battery Replacement</td><td></td></tr><tr><td colspan="2">☐ Dry Gas Regulator Replacement</td><td></td></tr><tr><td colspan="2">☐ Breath Tube Replacement</td><td></td></tr><tr><td colspan="2">☐ Other _____</td><td></td></tr></tbody></table>			Maintenance By PN		Date <u>5/29/2025</u>	☒ Battery Replacement			☐ Dry Gas Regulator Replacement			☐ Breath Tube Replacement			☐ Other _____																																					
Simulator	Serial #	Lot #/Exp																																																																							
0.050	MP5088	202406K 06/19/2026																																																																							
0.080	MP5089	202406L 06/19/2026																																																																							
0.200	MP5090	202406N 06/20/2026																																																																							
0.080 DGS	N/A	AG429602 10/22/2026																																																																							
Maintenance By PN		Date <u>5/29/2025</u>																																																																							
☒ Battery Replacement																																																																									
☐ Dry Gas Regulator Replacement																																																																									
☐ Breath Tube Replacement																																																																									
☐ Other _____																																																																									
Calibration Adjustment By _____			Department Inspection By <u>PN</u>																																																																						
Barometric Pressure Gauge _____ ID # _____ <table border="1"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #</th><th>Expiration</th></tr></thead><tbody><tr><td>0.000</td><td></td><td>N/A</td><td>N/A</td></tr><tr><td>0.040</td><td></td><td></td><td></td></tr><tr><td>0.100</td><td></td><td></td><td></td></tr><tr><td>0.200</td><td></td><td></td><td></td></tr><tr><td>0.300</td><td></td><td></td><td></td></tr><tr><td>0.080 DGS</td><td>N/A</td><td></td><td></td></tr></tbody></table> ☐ Post Calibration Adjustment Stability Checks <table border="1"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #</th><th>Expiration</th></tr></thead><tbody><tr><td>0.050</td><td></td><td></td><td></td></tr><tr><td>0.080</td><td></td><td></td><td></td></tr><tr><td>0.200</td><td></td><td></td><td></td></tr><tr><td>0.080 DGS</td><td>N/A</td><td></td><td></td></tr></tbody></table>			Simulator	Serial #	Lot #	Expiration	0.000		N/A	N/A	0.040				0.100				0.200				0.300				0.080 DGS	N/A			Simulator	Serial #	Lot #	Expiration	0.050				0.080				0.200				0.080 DGS	N/A			Barometric Pressure ID# <u>28427</u> Gauge <u>1010</u> Instrument <u>1007</u> Mouth Alcohol Solution Lot # <u>2025-A</u> Acetone Stock Solution Lot # <u>2024-B</u> <table border="1"><thead><tr><th>Simulator</th><th>Serial Number</th></tr></thead><tbody><tr><td>0.000</td><td>MP5086</td></tr><tr><td>Interferent</td><td>MP5087</td></tr><tr><td>0.050</td><td>MP5088</td></tr><tr><td>0.080</td><td>MP5089</td></tr><tr><td>0.200</td><td>MP5090</td></tr></tbody></table> Attachments <table border="1"><tbody><tr><td>☒ Form 41</td><td>☐ Post-Stability Checks</td></tr><tr><td>☒ Stability Checks</td><td>☐ Flow Calibration</td></tr><tr><td>☒ Calibration Certificate</td><td>☐ Form 40</td></tr><tr><td>☐ Calibration Adjustment</td><td>☐ Other _____</td></tr></tbody></table> ☒ Instrument Complies with Chapter 11D-8, FAC ☐ Instrument Does Not Comply with Chapter 11D-8, FAC ☒ Return to/Place into Evidentiary Use ☐ Remain Out of Evidentiary Use ☒ Conduct an Agency Inspection Before Evidentiary Use Digitally signed by Shayla Platt Taylor Gutschow Date: 2025.06.03 08:23:16 -04'00' Shayla Platt Date: 2025.06.03 09:20:07 -04'00' Tech Review / Date Admin Review / Date			Simulator	Serial Number	0.000	MP5086	Interferent	MP5087	0.050	MP5088	0.080	MP5089	0.200	MP5090	☒ Form 41	☐ Post-Stability Checks	☒ Stability Checks	☐ Flow Calibration	☒ Calibration Certificate	☐ Form 40	☐ Calibration Adjustment	☐ Other _____
Simulator	Serial #	Lot #	Expiration																																																																						
0.000		N/A	N/A																																																																						
0.040																																																																									
0.100																																																																									
0.200																																																																									
0.300																																																																									
0.080 DGS	N/A																																																																								
Simulator	Serial #	Lot #	Expiration																																																																						
0.050																																																																									
0.080																																																																									
0.200																																																																									
0.080 DGS	N/A																																																																								
Simulator	Serial Number																																																																								
0.000	MP5086																																																																								
Interferent	MP5087																																																																								
0.050	MP5088																																																																								
0.080	MP5089																																																																								
0.200	MP5090																																																																								
☒ Form 41	☐ Post-Stability Checks																																																																								
☒ Stability Checks	☐ Flow Calibration																																																																								
☒ Calibration Certificate	☐ Form 40																																																																								
☐ Calibration Adjustment	☐ Other _____																																																																								
Notes/Suggested Service: _____ _____ _____ _____ _____ _____																																																																									

STABILITY CHECKS

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000
05/30/2025
Software: 8100.27
SN 80-000779

Test	g/210L	Time
Air Blank	0.000	10:47
Control Test	0.052	10:47
Air Blank	0.000	10:48
Control Test	0.053	10:49
Air Blank	0.000	10:49
Control Test	0.052	10:50
Air Blank	0.000	10:50
Control Test Stats		
Average	0.0523	
Std Dev	0.0006	
Rel Std Dev(%)	1.1032	

Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000
05/30/2025
Software: 8100.27
SN 80-000779

Test	g/210L	Time
Air Blank	0.000	10:53
Control Test	0.081	10:53
Air Blank	0.000	10:54
Control Test	0.080	10:55
Air Blank	0.000	10:55
Control Test	0.081	10:56
Air Blank	0.000	10:56
Control Test Stats		
Average	0.0817	
Std Dev	0.0006	
Rel Std Dev(%)	0.7157	

Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000
05/30/2025
Software: 8100.27
SN 80-000779

Test	g/210L	Time
Air Blank	0.000	11:00
Control Test	0.200	11:00
Air Blank	0.000	11:01
Control Test	0.200	11:02
Air Blank	0.000	11:02
Control Test	0.201	11:03
Air Blank	0.000	11:03
Control Test Stats		
Average	0.2003	
Std Dev	0.0006	
Rel Std Dev(%)	0.2882	

Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000
05/30/2025
Software: 8100.27
SN 80-000779

Test	g/210L	Time
Air Blank	0.000	11:06
Control Test	0.080	11:06
Air Blank	0.000	11:06
Control Test	0.079	11:07
Air Blank	0.000	11:07
Control Test	0.079	11:08
Air Blank	0.000	11:08
Control Test Stats		
Average	0.0793	
Std Dev	0.0006	
Rel Std Dev(%)	0.7277	

DGS

Operator's Signature

Florida Department of Law Enforcement Alcohol Testing Program

DEPARTMENT INSPECTION REPORT - INTOXILYZER 8000

Agency: MADISON COUNTY SO
Time of Inspection: 13:22

Date of Inspection: 05/30/2025

Serial Number: 80-000779
Software: 8100.27

Check or Test	YES	NO	Check or Test	YES	NO
Diagnostic Check (Pre-Inspection): OK	Yes		Date and/or Time Adjusted		No
Minimum Sample Volume Check: OK	Yes		Barometric Pressure Sensor Check: OK	Yes	
Alcohol Free Subject Test: 0.000	Yes		Mouth Alcohol Test: Slope Not Met	Yes	
Interferent Detect Test: Interferent Detect	Yes		Diagnostic Check (Post-Inspection): OK	Yes	

Alcohol Free Test (g/210L)	0.05g/210L Test (g/210L) Lot#:202406K Exp: 06/19/2026	0.08g/210L Test (g/210L) Lot#:202406L Exp: 06/19/2026	0.20g/210L Test (g/210L) Lot#:202406N Exp: 06/20/2026	0.08 g/210L Dry Gas Std Test (g/210L) Lot#:AG429602 Exp: 10/22/2026
/ 0.000	0.052	0.081	0.194	0.080
/ 0.000	0.052	0.079	0.194	0.079
/ 0.000	0.052	0.078	0.195	0.080
/ 0.000	0.051	0.078	0.195	0.080
/ 0.000	0.052	0.078	0.194	0.080
/ 0.000	0.052	0.078	0.194	0.080
/ 0.000	0.053	0.078	0.194	0.080
/ 0.000	0.052	0.078	0.194	0.080
/ 0.000	0.053	0.078	0.194	0.080
/ 0.000	0.053	0.079	0.194	0.079

Standard Deviations	0.0006	0.0009	0.0004	0.0004
---------------------	--------	--------	--------	--------

Average Standard Deviation of 0.05, 0.08 and 0.20 g/210L Tests: 0.0005 Number of Simulators Used: 5

Remarks:

00: Ambient Fail: TURNED ON FAN AND CLEARED AREA.


. Int Det: Ambient Fail:

*Inspector in training used highly scented Hand Lotion in room.
Turned two fans on to ensure room was cleared of
Scent. - PN 5/30/2025*

Repeated the Alcohol-Free & Interferent Detect test. Each had an ambient fail on the first attempt. Corrective action (see above) resulted in passing test. PN 5/30/2025


The above instrument complies (☒) does not comply (☐) with Chapter 11D-8, FAC.

I certify that I performed this inspection in accordance with the provisions of Chapter 11D-8, FAC.

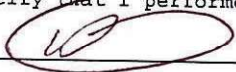

PHIL NICODEMO
Signature and Printed Name

05/30/2025
Date

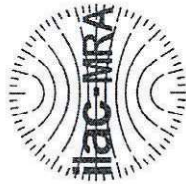

Calibration Certificate

Florida Department of Law Enforcement
Alcohol Testing Program
2331 Phillips Road
Tallahassee, FL 32308

This is to certify the calibration of Intoxilyzer 8000 serial number 80-000779, manufactured by CMI, Inc. was calibrated in accordance with FDLE/ATP Form 36 - Department Inspection Procedures - Intoxilyzer 8000.

Serial Number:	<u>80-000779</u>	UNCERTAINTY* $\pm$
Owning Agency:	<u>MADISON COUNTY SO</u>	0.050 g/ 210 L 0.004
Calibration Date:	<u>05/30/2025</u>	0.080 g/ 210 L 0.004
Calibration Time:	<u>13:22</u>	0.200 g/ 210 L 0.007
		0.080 g/ 210 L Dry Gas Control 0.005

All results are reported in g/ 210 L.

Bias is limited by calibration acceptance criteria. All calibration results must be within ± 0.005 or 5%, whichever is greater, of the target alcohol concentration.

*Uncertainty is based on fleet-wide data and is expressed to a 99.73% level of confidence ($k=3$).

The instrument results before and after any adjustment are found in the associated pre and post stability checks.

TRACEABILITY INFORMATION

This instrument was calibrated using solutions prepared by Alcohol Countermeasure Systems, Inc. (ACS). ACS prepared and certified these CRMs in accordance with ISO 17034 and ISO/ IEC 17025 Standards.

Simulator temperatures are traceable to NIST. Simulator temperatures are checked with NIST traceable digital thermometers calibrated by Precision Metrology in accordance with ISO/ IEC 17025 standards.

Dry gas control measurements are traceable to NIST through the use of CRMs supplied by an accredited CRM supplier. The supplier of dry gas standard controls prepared and certified the CRMs in accordance with ISO Guide 34 and ISO/ IEC 17025 standards.

This document shall not be reproduced except in full, without written approval of the Florida Department of Law Enforcement Alcohol Testing Program.

05/30/2025

Date


PHIL NICODEMO,
Department Inspector

FDLE/ATP Form 69 October 2024
Issuing Authority: Alcohol Testing Program

Service • Integrity • Respect • Quality



INSTRUMENT PROCESSING SHEET

Agency Madison County SOS/N 80-000779Florida Department of
Law EnforcementDate In 12/30/2024 DI Completion Date 01/08/2025☐ Ship ☒ P/U ☐ H/D ☐ CMI ☐ EE

Intake By <u>ALL</u> Date <u>12/30/2024</u>		Quality Checks By <u>DA</u> Date <u>01/08/2025</u>		Flow Calibration By _____ Date _____																																																													
☒ Annual ☐ Registration ☐ Return from CMI / EE Visual Inspection: ☒ Case ☐ Handle ☒ Keyboard ☐ Dry Gas Shelf ☒ Feet ☐ Breath Tube ☒ Ports ☐ Screws Tight Other Equipment/ Accessories: ☐ Power cord ☐ Printer Cable ☐ Static Bag ☐ 12V DC Cable Notes: <u>dropoff, no box</u>		☒ Breath Tube Screen ☒ Replace External O-Rings ☒ Instrument Set Up Verified ☒ R-Value <u>226</u> ☒ Flow Verification (L/s) Flow Column # <u>ATP102</u> 32 mm <u>0.160</u> (.139 - .169) 36 mm <u>0.175</u> (.156 - .190) 53 mm <u>0.246</u> (.228 - .278) 103 mm <u>0.523</u> (.447 - .547) ☒ Barometric Pressure Check Gauge ID # <u>30793</u> ☒ Stability Checks		☐ Flow Column # _____ ☐ 5L/min – 17mm ☐ 15L/min – 53mm ☐ 30L/min – 103mm ☐ R-Value _____ ☐ Post Calibration Verification (L/s) Flow Column # _____ 32 mm _____ (.139 - .169) 36 mm _____ (.156 - .190) 53 mm _____ (.228 - .278) 103 mm _____ (.447 - .547)																																																													
		<table border="1" style="width:100%"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #/Exp</th></tr></thead><tbody><tr><td>0.050</td><td>MP5088</td><td>202406K 06/19/2026</td></tr><tr><td>0.080</td><td>MP5089</td><td>202406L 06/19/2026</td></tr><tr><td>0.200</td><td>MP5090</td><td>202406N 06/20/2026</td></tr><tr><td>0.080 DGS</td><td>N/A</td><td>AG429602 10/22/2026</td></tr></tbody></table>		Simulator	Serial #	Lot #/Exp	0.050	MP5088	202406K 06/19/2026	0.080	MP5089	202406L 06/19/2026	0.200	MP5090	202406N 06/20/2026	0.080 DGS	N/A	AG429602 10/22/2026	Maintenance By _____ Date _____ ☐ Battery Replacement ☐ Dry Gas Regulator Replacement ☐ Breath Tube Replacement ☐ Other _____																																														
Simulator	Serial #	Lot #/Exp																																																															
0.050	MP5088	202406K 06/19/2026																																																															
0.080	MP5089	202406L 06/19/2026																																																															
0.200	MP5090	202406N 06/20/2026																																																															
0.080 DGS	N/A	AG429602 10/22/2026																																																															
Calibration Adjustment By _____				Department Inspection By <u>DA</u>																																																													
Barometric Pressure Gauge _____ ID # _____ <table border="1" style="width:100%"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #</th><th>Expiration</th></tr></thead><tbody><tr><td>0.000</td><td></td><td>N/A</td><td>N/A</td></tr><tr><td>0.040</td><td></td><td></td><td></td></tr><tr><td>0.100</td><td></td><td></td><td></td></tr><tr><td>0.200</td><td></td><td></td><td></td></tr><tr><td>0.300</td><td></td><td></td><td></td></tr><tr><td>0.080 DGS</td><td>N/A</td><td></td><td></td></tr></tbody></table> ☐ Post Calibration Adjustment Stability Checks <table border="1" style="width:100%"><thead><tr><th>Simulator</th><th>Serial #</th><th>Lot #</th><th>Expiration</th></tr></thead><tbody><tr><td>0.050</td><td></td><td></td><td></td></tr><tr><td>0.080</td><td></td><td></td><td></td></tr><tr><td>0.200</td><td></td><td></td><td></td></tr><tr><td>0.080 DGS</td><td>N/A</td><td></td><td></td></tr></tbody></table>				Simulator	Serial #	Lot #	Expiration	0.000		N/A	N/A	0.040				0.100				0.200				0.300				0.080 DGS	N/A			Simulator	Serial #	Lot #	Expiration	0.050				0.080				0.200				0.080 DGS	N/A			Barometric Pressure ID# <u>30793</u> Gauge <u>1021</u> Instrument <u>1020</u> Mouth Alcohol Solution Lot # <u>2024-A</u> Acetone Stock Solution Lot # <u>2023-B</u> <table border="1" style="width:100%"><thead><tr><th>Simulator</th><th>Serial Number</th></tr></thead><tbody><tr><td>0.000</td><td>MP5086</td></tr><tr><td>Interferent</td><td>MP5087</td></tr><tr><td>0.050</td><td>MP5088</td></tr><tr><td>0.080</td><td>MP5089</td></tr><tr><td>0.200</td><td>MP5090</td></tr></tbody></table>		Simulator	Serial Number	0.000	MP5086	Interferent	MP5087	0.050	MP5088	0.080	MP5089	0.200	MP5090
Simulator	Serial #	Lot #	Expiration																																																														
0.000		N/A	N/A																																																														
0.040																																																																	
0.100																																																																	
0.200																																																																	
0.300																																																																	
0.080 DGS	N/A																																																																
Simulator	Serial #	Lot #	Expiration																																																														
0.050																																																																	
0.080																																																																	
0.200																																																																	
0.080 DGS	N/A																																																																
Simulator	Serial Number																																																																
0.000	MP5086																																																																
Interferent	MP5087																																																																
0.050	MP5088																																																																
0.080	MP5089																																																																
0.200	MP5090																																																																
Notes/Suggested Service: _____ _____ _____ _____ _____ _____				Attachments ☒ Form 41 ☐ Post-Stability Checks ☒ Stability Checks ☐ Flow Calibration ☒ Calibration Certificate ☐ Form 40 ☐ Calibration Adjustment ☐ Other _____																																																													
				☒ Instrument Complies with Chapter 11D-8, FAC ☐ Instrument Does Not Comply with Chapter 11D-8, FAC ☒ Return to/Place into Evidentiary Use ☐ Remain Out of Evidentiary Use ☒ Conduct an Agency Inspection Before Evidentiary Use Taylor Gutschow Digitally signed by Taylor Gutschow Date: 2025.01.14 14:53:39 -05'00' Shayla Platt Digitally signed by Shayla Platt Date: 2025.01.14 15:13:38 -05'00' Tech Review / Date Admin Review / Date																																																													

Stability Checks 80-000779 1/8/25 DA

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000779
01/08/2025
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	12:24
Control Test	0.052	12:24
Air Blank	0.000	12:25
Control Test	0.053	12:25
Air Blank	0.000	12:26
Control Test	0.053	12:27
Air Blank	0.000	12:27
Control Test Stats		
Average	0.0527	
Std Dev	0.0006	
Rel Std Dev (%)	1.0962	



Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000779
01/08/2025
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	12:28
Control Test	0.082	12:29
Air Blank	0.000	12:30
Control Test	0.082	12:30
Air Blank	0.000	12:31
Control Test	0.082	12:32
Air Blank	0.000	12:32
Control Test Stats		
Average	0.0820	
Std Dev	0.0000	
Rel Std Dev(%)	0.0000	


wet



Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000779
01/08/2025
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	12:34
Control Test	0.202	12:34
Air Blank	0.000	12:35
Control Test	0.202	12:35
Air Blank	0.000	12:36
Control Test	0.202	12:37
Air Blank	0.000	12:37
Control Test Stats		
Average	0.2020	
Std Dev	0.0000	
Rel Std Dev(%)	0.0000	

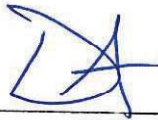


Operator's Signature

MADISON COUNTY SO
Intoxilyzer - Alcohol Analyzer
Model 8000 SN 80-000779
01/08/2025
Software: 8100.27

Test	g/210L	Time
Air Blank	0.000	12:39
Control Test	0.081	12:39
Air Blank	0.000	12:40
Control Test	0.082	12:40
Air Blank	0.000	12:41
Control Test	0.081	12:41
Air Blank	0.000	12:41
Control Test Stats		
Average	0.0813	
Std Dev	0.0006	
Rel Std Dev(%)	0.7099	

DGS

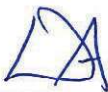

Operator's Signature

Florida Department of Law Enforcement Alcohol Testing Program

DEPARTMENT INSPECTION REPORT - INTOXILYZER 8000

Agency: MADISON COUNTY SO
Time of Inspection: 14:22

Date of Inspection: 01/08/2025

Serial Number: 80-000779
Software: 8100.27

Check or Test	YES	NO	Check or Test	YES	NO
Diagnostic Check (Pre-Inspection): OK	Yes		Date and/or Time Adjusted	Yes	
Minimum Sample Volume Check: OK	Yes		Barometric Pressure Sensor Check: OK	Yes	
Alcohol Free Subject Test: 0.000	Yes		Mouth Alcohol Test: Slope Not Met	Yes	
Interferent Detect Test: Interferent Detect	Yes		Diagnostic Check (Post-Inspection): OK	Yes	

Alcohol Free Test (g/210L)	0.05g/210L Test (g/210L) Lot#:202406K Exp: 06/19/2026	0.08g/210L Test (g/210L) Lot#:202406L Exp: 06/19/2026	0.20g/210L Test (g/210L) Lot#:202406N Exp: 06/20/2026	0.08 g/210L Dry Gas Std Test (g/210L) Lot#:AG429602 Exp: 10/22/2026
0.000	0.051	0.080	0.199	0.082
0.000	0.052	0.080	0.199	0.081
0.000	0.052	0.080	0.198	0.081
0.000	0.052	0.080	0.201	0.081
0.000	0.051	0.081	0.201	0.081
0.000	0.051	0.080	0.199	0.081
0.000	0.051	0.081	0.199	0.081
0.000	0.051	0.081	0.200	0.081
0.000	0.051	0.081	0.200	0.081
0.000	0.051	0.081	0.201	0.081

Standard Deviations	0.0004	0.0005	0.0010	0.0003
---------------------	--------	--------	--------	--------


Average Standard Deviation of 0.05, 0.08 and 0.20 g/210L Tests: 0.0005 Number of Simulators Used: 5

Remarks:

Time-Date changed.TIME/DATE ACCIDENTALLY TYPED IN

The above instrument complies (☒) does not comply (☐) with Chapter 11D-8, FAC.

I certify that I performed this inspection in accordance with the provisions of Chapter 11D-8, FAC.

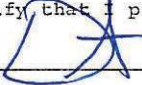
Signature and Printed Name

DESTINEE N ARMSTRONG

01/08/2025
Date



Calibration Certificate

Florida Department of Law Enforcement
Alcohol Testing Program
2331 Phillips Road
Tallahassee, FL 32308

This is to certify the calibration of Intoxilyzer 8000 serial number 80-000779, manufactured by CMI, Inc. was calibrated in accordance with FDLE/ATP Form 36 - Department Inspection Procedures - Intoxilyzer 8000.

Serial Number:	<u>80-000779</u>	UNCERTAINTY* $\pm$
Owning Agency:	<u>MADISON COUNTY SO</u>	0.050 g/ 210 L 0.004
Calibration Date:	<u>01/08/2025</u>	0.080 g/ 210 L 0.004
Calibration Time:	<u>14:22</u>	0.200 g/ 210 L 0.007
		0.080 g/ 210 L Dry Gas Control 0.005

All results are reported in g/ 210 L.

Bias is limited by calibration acceptance criteria. All calibration results must be within ± 0.005 or 5%, whichever is greater, of the target alcohol concentration.

*Uncertainty is based on fleet-wide data and is expressed to a 99.73% level of confidence ($k=3$).

The instrument results before and after any adjustment are found in the associated pre and post stability checks.

TRACEABILITY INFORMATION

This instrument was calibrated using solutions prepared by Alcohol Countermeasure Systems, Inc. (ACS). ACS prepared and certified these CRMs in accordance with ISO 17034 and ISO/ IEC 17025 Standards.

Simulator temperatures are traceable to NIST. Simulator temperatures are checked with NIST traceable digital thermometers calibrated by Precision Metrology in accordance with ISO/ IEC 17025 standards.

Dry gas control measurements are traceable to NIST through the use of CRMs supplied by an accredited CRM supplier. The supplier of dry gas standard controls prepared and certified the CRMs in accordance with ISO Guide 34 and ISO/ IEC 17025 standards.

This document shall not be reproduced except in full, without written approval of the Florida Department of Law Enforcement Alcohol Testing Program.

Destinee
Armstrong
Digitally signed by Destinee
Armstrong
Date: 2025.01.14 08:59:48
+05'00'

01/08/2025

Date

DESTINEE N ARMSTRONG,
Department Inspector

FDLE/ATP Form 69 October 2024

Issuing Authority: Alcohol Testing Program

Service • Integrity • Respect • Quality